

Sample Preparation Guide FS




Experimental Design: The FlowSight system can quantify the intensity, specific location, and distribution of signals within tens of thousands of cells per sample. The system can perform most flow cytometric assays, but the best applications take advantage of the technology's imaging capabilities to discriminate subtle morphological changes within individual cells and cell populations.

FlowSight


- Choice of Cell Type:** The particle size should be less than 60um using 20x magnification.

Basic example (2um/pixel)

THP-1 (~20 um diameter)

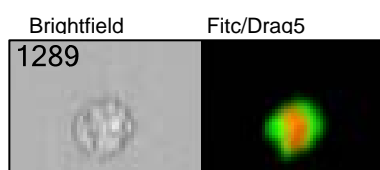


Blood Cell (~8um diameter)

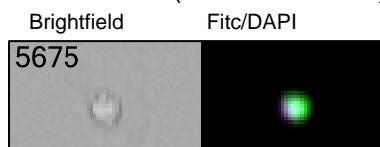


QI example (1um/pixel)

THP-1

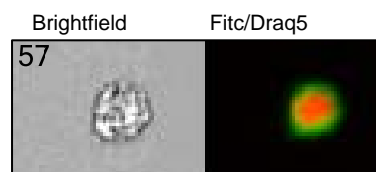

Blood Cell

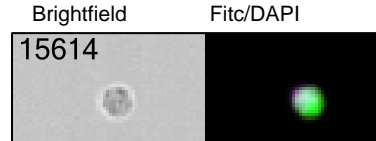

- Final Sample Concentration and Volume:** At least 1 million cells in 50 μL (2×10^7 cells/ml) in PBS/2%FBS in a 1.5mL siliconized microcentrifuge tube.
- Protocols:** In general, any established labeling protocol used for flow cytometry will work with the FlowSight (see *Current Protocols in Cytometry* for general labeling techniques). Stain cells on ice in the presence of azide when possible to reduce non-specific capping of antibody. Use polypropylene tubes, preferably siliconized, to process samples.
- Choice of Fluorochromes:** Choose fluorochromes that are excited by the lasers in your FlowSight (405,488,642nm are most common). Use the chart on p.3 or look online for a spectra viewer that will help you plan which dyes will work the best.
- Compensation:** Have a sample of cells each labeled with a single-color for each fluorochrome used (i.e. FITC only cells, PE only cells, etc.).
- Cell Aggregation:** Minimize aggregation problems by straining the sample through a 70um nylon mesh strainer, or by using an anti-clumping buffer such as EDTA or Accumax prior to fixation.
- Fixation:** If fixation is desired, thoroughly fix cells with 1% PFA on ice for 20 min.
- Number of samples:** No more than 30 total for feasibility experiments. Please limit the samples to the following; Positive and Negative **biologic controls**, **compensation controls**, and **experiment samples**.

Sample Preparation Guide



9. Brightness of Stain and Stain Balancing: Quantifying the location and distribution of signals in an image is a demanding task that requires optimized labeling. Below are a few suggestions to help design the experiment:

- Try to achieve at least a full log shift in fluorescence, as measured by FACS.
- Use the brightest dye for the antigen with the smallest copy number.
- The brightness of probes can be independently controlled by changing the laser power. However, data quality is enhanced when the brightness levels of all probes excited off a single laser are balanced to within a log of each other. Probe balancing avoids the saturation of bright stains when they are combined with dim stains in the same sample.

10. Shipping recommendations:

- Include the [data acquisition form](#) and data that verifies the experiment worked.
- Samples must be fixed and non-pathogenic.
- Wrap sample tubes in Parafilm.
- Insulate by placing in 15ml conical rack with Styrofoam lid. Pack in Styrofoam lined shipping box with refrigerant packs and paper. For winter delivery, use room temperature packs to prevent freezing of sample. For summer delivery use frozen packs.
- Label outside of box 'Do Not Freeze'.
- Email the package tracking number to your Amnis contact person.

11. International Shipments

To ship non-viable (fixed) material internationally to Amnis, a written statement for US Customs and Border Protection, Department of Homeland Security **must** accompany the shipment. ***This written statement must:***

- Be addressed to; US Customs and Border Protection, Department of Homeland Security.
- Be an original copy on institutional letterhead signed by the laboratory worker responsible for preparing the samples.
- Identify the material and name of the species from which the material was derived.
- state that the animals from which the material was derived:
 - a. have not been exposed to, or inoculated with, any livestock or poultry disease agents exotic to the United States, and
 - b. did not originate from a facility where work with exotic disease agents affecting livestock or avian species is conducted
- State that the material is non-viable
- Be placed in an envelope addressed to 'US Customs and Border Protection, Department of Homeland Security' and attached to the **outside** of the shipping box.

Shipping Address:

Amnis Corporation

645 Elliott Avenue West, Suite 100

Seattle WA 98119

Pn: (855) 240-8045

Fx: (206) 576-6895

Email : applicationsupport@amnis.com
www.amnis.com





Sample Preparation Guide

FS Quick Start Guide: Sample Prep Guide

		Excitation Laser (nm)							
Ch	Band (nm)	405	488	561	642	785	Used	Ch	
1	435-505 (457/45)	BRIGHTFIELD							1
2	505-560 (532/55)		FITC, AF488, GFP, YFP, DyLight488, PKH67, Syto13, SpectrumGreen, LysoTrackerGreen, MitoTrackerGreen					2	
3	560-595 (577/35)		PE*, PKH26, Cy3, DSRed, CellMask/CellTracker/ SYTOX Orange	PE*, AF546, Cy3*, DyLight550, PKH26, DSRed, SpectrumOrange, MitoTrackerOrange				3	
4	595-642 (610/30)		PE-TexRed*, ECD*, PE- AF610*, PI*, RFP, eFluor625*	AF568*, AF594*, AF610*, Cy3*, DyLight594*, PE- TexRed*, ECD*, TexRed*, PE-AF610*, RFP, mCherry*, PI*				4	
5	642-745 (702/86)		PE-Cy5*, PE-AF647*, PerCP*, PerCP- Cy5.5*, 7AAD* eFluor650* FuraRedlo	PE-Cy5*, PE-AF647* 7AAD*				5	
6	745-800 (772/55)		PE-Cy7*, PE-AF750*, Draq5*, LDS751*	PE-Cy7*, PE-AF750*, Draq5*, LDS751*		SSC		6	
7	435-505 (457/45)	DAPI, BV421, AF405, Hoechst, PacBlue, CascadeBlue, eFluor450, DyLight405, CFP, LIVE/DEAD Violet						7	
8	505-560 (532/55)	BV510, PacOrange, CascadeYellow, AF430, eFluor525, QD525,						8	
9	560-595 (577/35)	BRIGHTFIELD							9
10	595-642 (610/30)	QD625*, eFluor625*, BV605						10	
11	642-745 (702/86)	QD705*, eFluor700, BV711			AF647, AF660, AF680, DRAQ5*, APC, Cy5, DyLight649, DyLight680, PE-AF647*, PE-Cy5*, PerCP*, PerCP-Cy5.5*			11	
12	745-800 (772/55)	QD800*, BV786			APC-Cy7, APC-AF750, APC-H7, APC-Cy7, Cy7, AF750, DyLight750, PE- Cy7*, PE-AF750*,			12	

Recommended dyes (based on optimal excitation and detection channels) are in boldface.

1 laser (488) system: Ideal dyes are AF488, PE, PE-TxRed, PE-Cy5, SSC-Ch6,

2 laser (488,642) system: Ideal dyes are AF488, PE, PE-TxRed, SSC-Ch6, and AF647, APC Cy7

3 laser (405,488,642) system: Ideal dyes are AF488, PE, PE-TxRed, SSC-Ch6, and DAPI, AF647, APC Cy7

Band-pass filter values assume 3laser configuration.



Part of EMD Millipore