

Department Of Veterans Affairs  
Decentralized Hospital Computer Program

# **LABORATORY VITEK LITERAL INTERFACE LA\*5.2\*12 PATCH DOCUMENTATION**

Version 5.2

April 1996

Information Resources Management Field Office  
Birmingham, Alabama



# Preface

The Decentralized Hospital Computer Program (DHCP) Laboratory Literal Vitek Interface LA\*5.2\*12 Patch Documentation is designed to provide the Department of Veterans Affairs Medical Center (DVAMC), Information Resources Management Service (IRM), and the Laboratory Microbiology Staff with the necessary technical information required to efficiently and effectively implement, maintain, and manage the bi-directionally literal interface and the Vitek Automated Instrument to DHCP through the Laboratory System Interface (LSI). The LA\*5.2\*12 Patch Documentation also provides sufficient information about DHCP downloading data to Vitek and Vitek uploading data to DHCP.



# Table of Contents

Introduction .....	7
New Features:.....	8
References: .....	8
Documentation Notations:.....	8
Keywords and Phrases: .....	9
Pre-Installation Information .....	11
Staffing Resources: .....	11
Test Account:.....	11
Routines: .....	11
Disk Space:.....	12
New Globals: .....	12
DHCP Software Requirements: .....	12
Installation Requirements:.....	12
Vitek Computer Requirements: .....	12
New Vitek Options:.....	13
New DHCP Files:.....	13
MICRO INSTRUMENT SET-UP file (#61.38) .....	13
LABORATORY INSTRUMENT CODE (LIC) file (#61.39) .....	14
LA*5.2*12 Patch Implementation Checklist .....	15
VITEK to DHCP Laboratory System Interface (LSI) Cable Wiring.....	17
bioMerieux Vitek Computer Setup .....	19
Ports Setup.....	19
Protocol Setup .....	20
LA*5.2*12 Patch Installation.....	21
Post-Installation Information .....	26
Laboratory Microbiology Staff.....	26
Print DHCP and Vitek Code Entries .....	26
Editing DHCP Files .....	28
LABORATORY INSTRUMENT CODE FILE file (#61.39) .....	28
LOAD/WORK LIST file (#68.2).....	31
AUTO INSTRUMENT file (#62.4) .....	32
MICRO INSTRUMENT SET-UP file (#61.38) .....	33
Vitek Literal Verification Process .....	36
Troubleshooting .....	42

## Table of Contents

# Introduction

Approximately 50 VAMC facilities have purchased the bioMerieux Vitek Analyzer to perform automated microbiology laboratory testing. The bioMerieux Vitek Analyzer performs bacterial identification and susceptibility profiles. The main benefit from having an interface from the analyzer to a Laboratory Information System (LIS) is the improvement in overall efficiency of information handling. Data processed on the bioMerieux Vitek and in the DHCP Laboratory Database can be transferred electronically between both computer systems quickly and more efficiently than manual transfers. Manual data entry is not only labor intensive, but prone to transcription errors.

The bioMerieux Vitek has been interfaced to the DHCP Laboratory software for several years in the uni-directional and bi-directional modes. Previously the data stream from the bioMerieux Vitek was encoded, that is, the data coded is in a non human readable format.

Within the last five years bioMerieux Vitek released a new version of software using a literal interface (Literal PRM). In the last couple of years bioMerieux has introduced an additional software update, the bioLiaison incorporating the Literal PRM Interface with Graphical User Interface or GUI Version. The interface that is included in this update can be used with both the bioMerieux Vitek Literal PRM mode and the bioLiaison GUI Version. The bioMerieux Vitek, Inc., has indicated that the newer updates will be for the Literal PRM Interface or GUI version, not the encoded interface. There will be some updating of the encoded interface but eventually will be phased out.

This project encompasses all necessary activities involved with the interfacing of the bioMerieux Vitek automated microbiology analyzer to DHCP Laboratory software. These activities include:

- ♦Transmission of DHCP data to the bioMerieux Vitek analyzer in a format which is acceptable to the Vitek
- ♦Accepting result data as it is transmitted from the bioMerieux Vitek and interpreting the result data
- ♦Presenting the result data to DHCP Laboratory software in a format suitable for user verification of tests results

## New Features:

The DHCP Laboratory Vitek Literal Interface includes a Translation Table file, LABORATORY INSTRUMENT CODE (LIC) file (#61.39). The LIC file (#61.39) replaces the entries in the card entries that are stored in the AUTO INSTRUMENT file (#62.4).

The MICRO INSTRUMENT SET-UP file (#61.38) provides the site greater flexibility for setting up wild cards fields for the download, and whether to enable the LABORATORY TEST file (#60) culture ID Prefix Field.

## References:

Bidirectional Computer Interface Specification for bioLiaison Vitek, Part Number 510703-1, bioMerieux Vitek, 595 Anglum Drive, Hazelwood, MO. 63042-2395, Phone: 314-731-8500, Fax: 314-731-8800.

Bidirectional Computer Interface Operator's Manual, 510710-1, REV 1094 bioMerieux Vitek, 595 Anglum Drive, Hazelwood, MO. 63042-2395, Phone: 314-731-8500, Fax: 314-731-8800.

DHCP Laboratory Technical Manual, V. 5.2, October 1994, pp. 114-143, Dept. of Veterans Affairs (DVA), Veterans Health Administration (VHA), Office of the Chief Information Officer (OCIO)

## Documentation Notations:

- ♦Computer dialogue appears in Courier font, no larger than 10 points.  
**Example:** (Courier font 10 points).
- ♦User response appears in **boldface type** Courier font, no larger than 10 points.
- ♦After users have completed entering information the <**RET**> symbol appears in boldface type.
- ♦The underline appears when a user's response is required.



## Keywords and Phrases:

**ASCII Control characters:** ENQ, ACK, NAK, and EOT.

**Bi-directional Computer Interface (BCI):** Facilitates the exchange of data between the computer in a bioMerieux Vitek instrument and the DHCP computer.

**bioLiaison (LSN):** Refers to the Vitek patient database software.

**Download:** The process by which information is transferred from the host computer to the automated instrument.

**Firmware:** Broadly, the system software permanently stored in a computer's read-only memory (ROM) or elsewhere in the computer's circuitry. Firmware cannot be modified by the user.

**LIS:** Laboratory Information System

**LSI:** Laboratory System Interface

**Session:** Consists of a connection between the Vitek and the host to exchange information. A session can take place in only one direction at a time, where the initiator is the master and the other is the slave. ENQ initiates while EOT ends the communication/message.

**Translation Table:** Programming that converts data from one system to another facilitating electronic communication.

**Upload:** The transfer of information from the automated instrument to the host computer.

**Vitek Literal Interface Software:** Facilitates the exchange of data between the computer in a bioMerieux Vitek instrument and the DHCP computer.



# Pre-Installation Information

This section of the documentation provides pre-installation information that is necessary to ensure a successful installation of the LA\*5.2\*12 patch and implementation of the DHCP Laboratory Vitek Literal Interface.

## Staffing Resources:

1. Personnel with VA FileMan capabilities are required for editing the following files:
  - a) ANTIMICROBIAL SUSCEPTIBILITY file (#62.06)
  - b) AUTO INSTRUMENT file (#62.4)
  - c) LOAD/WORK LIST file (#68.2)
  - d) ETIOLOGY FIELD file (#61.2)
  - e) LABORATORY INSTRUMENT CODE file (#61.39)
  - f) MICRO INSTRUMENT SET-UP file (#61.38)
2. Personnel required for wiring Vitek cable to the LSI.
3. Personnel required to setup port and protocol on the Vitek.

## Test Account:

A test account is not applicable due to the data required for testing. It is highly recommended to set up a Test Patient in your live account for data verification.

## Routines:

LAMIVTL0 (Verification)	LAMIVTLC (LAH Builder)
LAMIVTL1 (Verification)	LAMIVTLP (LA Parser)
LAMIVTL2 (Verification)	LAMIVTLX (Handshake)
LAMIVTL3 (Verification)	LAMIVTLG (New LAGEN)
LAMIVTL4 (Verification)	LAMIVTLD (Download Routine)
LAMIVTL5 (Verification)	LAMIVTLB (Pre-Inits)
LAMIVTL6 (Verification)	LAMIVTLE (Post-Inits)

## Disk Space:

Minimal disk space growth will occur from the installation of the LA\*5.2\*12 patch.

## New Globals:

- ♦MICRO INSTRUMENT SET-UP file (#61.38) is ^LAB(61.38).
- ♦LABORATORY INSTRUMENT CODE file (#61.39) global is ^LAB(61.39).

## DHCP Software Requirements:

DHCP Software	Laboratory V. 5.2 Files
Laboratory V. 5.2	ANTIMICROBIAL SUSCEPTIBILITY file (#62.06)
VA FileMan V. 21	AUTO INSTRUMENT file (#62.4)
Kernel V. 8.0	ETIOLOGY FIELD file (#61.2)
	LABORATORY INSTRUMENT CODE file (#61.39)
	LOAD/WORK LIST file (#68.2)
	MICRO INSTRUMENT SET-UP file (#61.38)

## Installation Requirements:

The LA\*5.2\*12 Patch installation is using the Kernel Installation and Distribution Systems (KIDS), a new module in Kernel Version 8.0.

## Vitek Computer Requirements:

### **VITEK VTK R02.0 (Graphical User Interface (GUI) version) or greater**

The Vitek text-based software programs have been converted and updated to the new “graphical” platform.

### **Bi-directional Computer Interface (BCI) Version R020 or greater**

The BCI facilitates the exchange of data between the computer in a bioMerieux Vitek instrument and the DHCP computer.

## New Vitek Options:

1. Vitek Literal [LA VITEK LITERAL] option is used to access verify results. This option is added to the Microbiology [LRMI] menu by the KIDS installation process.
2. Antibiotic Code Print [LA PRINT ANTIBIOTIC] option is used to print the antibiotic code entries from the new LIC file (#61.39). This option is added to Microbiology Print [MICROBIOLOGY PRINT] MENU by the KIDS installation process.
3. Organism Code Print [LA PRINT ORGANISM] option is used to print the organism code entries from the new LIC file (#61.39). This option is added to Microbiology Print [MICROBIOLOGY PRINT] MENU by the KIDS installation process.

## New DHCP Files:

### MICRO INSTRUMENT SET-UP file (#61.38)

The MICRO INSTRUMENT SET-UP file (#61.38) is installed by KIDS. This file provides another dimension for performing tests on the Vitek. The Vitek only permits one test/accession number. This can be circumvented by entering a number between 0-9 in the prefix field of the LABORATORY TEST file (#60). Employing the logic that the Vitek will only accept up to a 6 digit accession number, and no accession shall exceed 99,999. The prefix number could be placed in front of the 5 digit accession. File (#61.38) also turns on the wild card fields for the Vitek.

**WARNING:** Use of these wild card fields may overwrite historical data in the Vitek.

**NOTE:** The development of the DHCP Laboratory Vitek Literal Interface project was done using an MSM (486) and ALPHA systems.

## LABORATORY INSTRUMENT CODE (LIC) file (#61.39)

The LABORATORY INSTRUMENT CODE (LIC) file (#61.39) (^LAB(61.39)) was created for the purpose of translating entries from information tables used by the laboratory automated instruments to entries located in DHCP files. For the DHCP Laboratory Vitek Literal Interface project, the LIC file (#61.39) facilitates the translation of the organism and antibiotic tables resident on the Vitek, to the DHCP Laboratory ETIOLOGY FIELD file (#61.2) and ANTIMICROBIAL SUSCEPTIBILITY file (#62.06) entries. The way this novel DHCP concept is accomplished is simple. The LIC file contains two multiple fields, the ORGANISM CODE field #1, subfile (#61.391) and the ANTIBIOTIC CODE field #2, subfile (#61.392). These are the main fields of the translator.

Using the organism as an example, once the Vitek identifies an organism it is uploaded across the interface to DHCP where it is matched to an entry in the ORGANISM CODE field #1 subfile (#61.391) of the LIC file (#61.39). The organism subfile also contains an interpretation entry, which describes the organism code and a local entry that points to the DHCP ETIOLOGY FIELD file (#61.2). The pointing action completes the translation loop between the Vitek and DHCP. The antibiotics tested against the organism by the Vitek are translated in the same manner as the organism.

The LABORATORY INSTRUMENT CODE file (#61.39) is installed by using KIDS. Some local editing of this file will be required. The LIC file (#61.39) comes with most of the codes needed to run the DHCP Vitek Literal Interface. However, additional fields have been added. The LIC file (#61.39) was edited by Barbara McCoy, Microbiology Supervisor at Muskogee VAMC. The organism and antibiotic codes from the Vitek were stored on a 3.5 floppy diskette. The codes were retrieved from the floppy diskette using the LAMIVTLC and LAMIVTLP upload routines. These two routines store the codes in an ^LA(9999) global. A utility routine was also created to enter the codes, via user approval, into the LIC file (#61.39).

# LA\*5.2\*12 Patch Implementation Checklist


This check list is provided to ensure a successful implementation of the DHCP Laboratory Vitek Literal Interface LA\*5.2\*12 patch. This checklist may be copied or removed from this section of the documentation for your convenience.

1. Connect cable from Vitek to LSI ☐
2. Set up bioMerieux Vitek Computer ports and protocol ☐
3. Run KIDS install ☐
4. Printout of antibiotic codes using the new  
Antibiotic Code Print [LA PRINT ANTIBIOTIC] option ☐
5. Printout of organism codes using the new  
Organism Code Print [LA PRINT ORGANISM] option ☐
6. Edit LIC file (#61.39) antibiotic field local entry pointing  
to the ANTIMICROBIAL SUSCEPTIBILITY file (#62.06) ☐
7. Edit LIC file (#61.39) organism field local entry pointing  
to the ETIOLOGY FIELD file (61.2) ☐
8. Edit LOAD/WORK LIST file (#68.2) ☐
9. Edit AUTO INSTRUMENT file (#62.4) ☐
10. Edit MICRO INSTRUMENT SET-UP file (#61.38) ☐






# VITEK to DHCP Laboratory System Interface (LSI) Cable Wiring

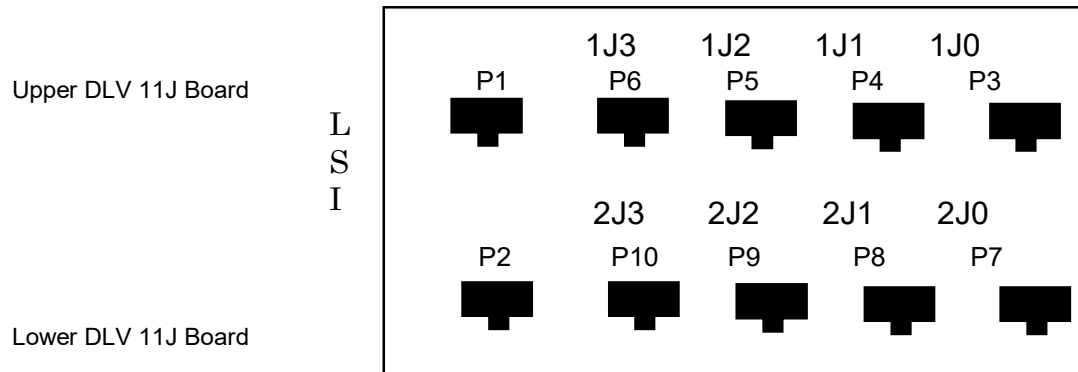

## LSI CABLES

All ports have been preset to 1200 baud by the factory. P1 is only to be used for the LSI-to -Host connection. The baud rates for ports P1 and P2 are hardwired to 1200 baud. All other ports may be site-configured from 150 to 9600 baud by wire wrapping or dip switch setting selected posts on the DLV11J board(s).

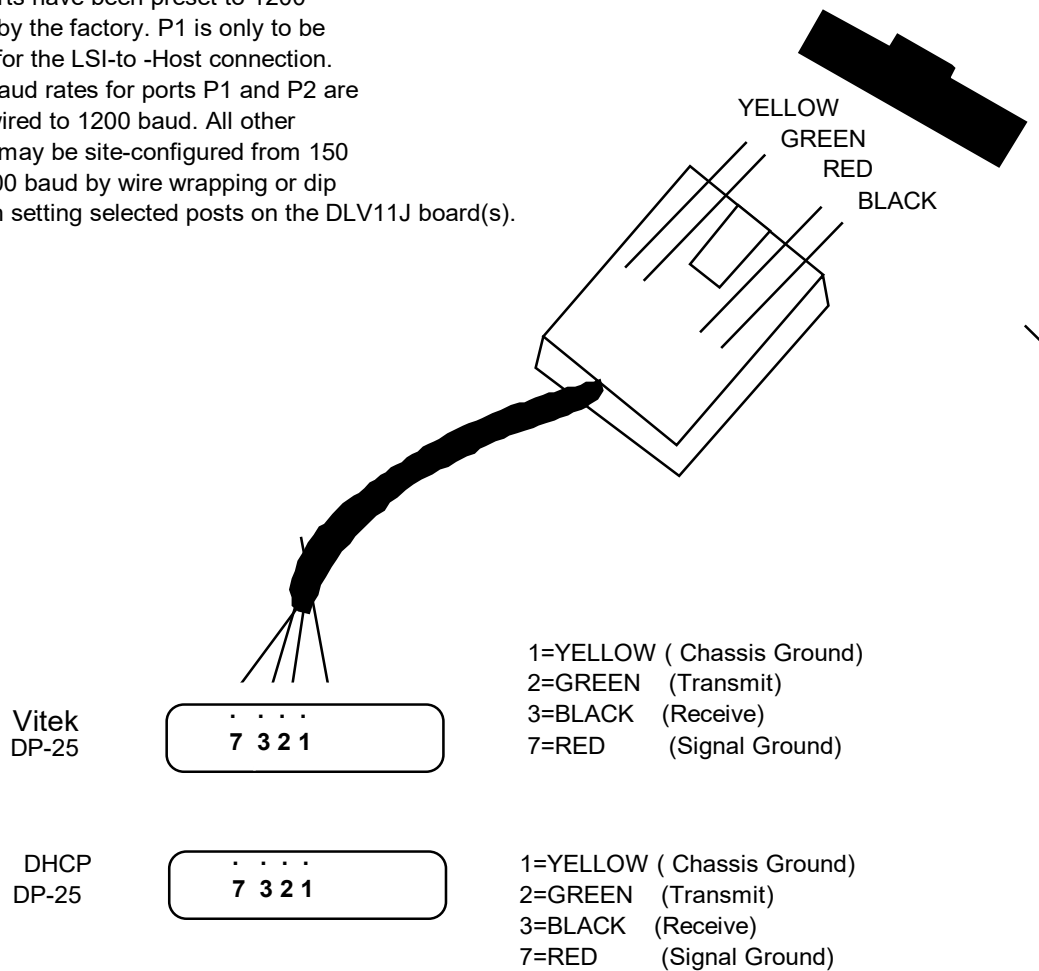



# bioMerieux Vitek Computer Setup

Use the following instructions for setting up the bioMerieux Vitek computer ports and protocol. The underline appears when a computer response is required.

## Ports Setup

1. From the bioMerieux screen:

Select - bioLiaison Main menu

2. From the bioLiaison Main menu:

Select - BCI

3. From the bioMerieux Vitek Bidirectional Computer Interface screen:

Select - PORTS

4. From the Single Select List screen:

Highlight - TTY0 and click

5. From the bioMerieuxx System Maintenance Screen; Change/Show Characteristics of a TTY, edit the following settings:

BAUD rate - Click on the LIST box and select 2400

PARITY - Click on the LIST box and select None

BITS per character - Click on the LIST box and select 8

Number of STOP BITS - Click on the LIST box and select 1

## Protocol Setup

1. From the bioLiaison Main menu:

Select - Protocol

2. From BCI Protocol Configuration (Port tty1) screen:

Select - Character Set ISO8859=1

Select - Field | (Select shift and bar key to enter the bar (|) symbol)

Select - Date (may leave blank)

Select - Time (may leave blank)

3. From the bioMerieuxx Literal Communication Protocol Parameter menu select the following parameter and enter setup data:

<CR><LF>		DELAYS	RETRIES	TIMEOUT
STX	DIS	Last Master <u>10</u>	<ENQ> Limit <u>3</u>	Checksum <u>10</u>
ETX	DIS	Inter Record <u>10</u>	<ENQ> Interval <u>10</u>	Host Response <u>10</u>
RS	DIS	Inter Message <u>2</u>	Checksum Limit <u>3</u>	
GS	<u>ENA</u>		Checksum Interval <u>10</u>	
ENQ	DIS			
EOT	DIS			

## LA\*5.2\*12 Patch Installation

The LA\*5.2\*12 Patch installation is using KIDS, a new module in Kernel V. 8.0. The LABORATORY INSTRUMENT CODE file (#61.39), MICRO INSTRUMENT SET-UP file (#61.38), and three new menu options are installed by this process. The LIC file (#61.39) comes with most of the entries needed to run the Vitek Literal Interface. However, some local editing of this will be required.

The LA\*5.2\*12 patch total installation time is less than 15 minutes.

Background jobs for Laboratory and other software packages are not affected by the LA\*5.2\*12 patch installation process.

All users may remain active on the system during the installation process.

### Example:

```
>D ^XUP<RET>
```

```
Setting up programmer environment
Terminal Type set to: C-VT100
```

```
Select OPTION NAME:      EVE          Systems Manager Menu<RET>
```

```
Select Systems Manager Menu Option:  Programmer Options <RET>
```

```
Select Programmer Options Option:    Kernel Installation & Distribution
System<RET>
```

```
Select Kernel Installation & Distribution System Option: MAILMan Menu<RET>
```

```
VA MailMan 7.1 service for HOAK.DAVID_R+@DEV.ISC-DALLAS.VA.GOV
You last used MailMan: 31 Mar 96 12:47
Your current banner is: Computer Programmer
```

```
Select MailMan Menu Option: READ a message<RET>
```

```
Read MAIL BASKET: IN//<RET>
```

```
LAST Message Number: 1050   Messages in BASKET: 1050<RET>
```

## LA\*5.2\*12 Installation Process

IN Basket Message: 1// 1050<RET>

Subj: LA\*5.2\*12 TEST 2 [#55472] 31 Mar 96 12:43 5506 Lines

From: HOAK,DAVID R. in 'IN' basket. Page 1

---

\$TXT Created by HOAK,DAVID R. at DEV.ISC-DALLAS.VA.GOV (DIFROM) on SUNDAY,  
03/31/96 at 12:43

\$END TXT

\$ROU XPDPINIT

XPDPINIT ;SFISC/RSD - Load a Packman Message using KIDS ; 6 Feb 95 14:48

; ;8.0;KERNEL;\*\*5\*\*;Jul 10, 1995

I '\$D(^XMB(3.9,+ \$G(XMZ),0)) W !!, "NO message to install!!" Q

I \$T(^XPDI PM)=" W !!, "KIDS doesn't exist!!" Q

G ^XPDI PM

\$END ROU XPDPINIT

\$TXT \$KIDS LA\*5.2\*12

\*\*INSTALL NAME\*\*

LA\*5.2\*12

"BLD",18,0)

LA\*5.2\*12^^0^2940927^n

"BLD",18,1,0)

Select MESSAGE Action: IGNORE (in IN basket)// X<RET>

Select PackMan function: **INSTALL/CHECK MESSAGE<RET>**

Warning: Installing this message will cause a permanent update of globals  
and routines and run the INIT. Do you really want to do this? NO// Y<RET>

Shall I preserve what is on disk in a separate back-up message ? YES//  
N<RET>

No backup message built.

Line 3 Message #55472 Unloading Routine ROU XPDPINIT  
LA\*5.2\*12

The following Entries already exist in the INSTALL file:

LA\*5.2\*12 Install Completed  
was loaded on Mar 31, 1996@12:48:58  
OK to continue? NO// YES<RET>.

Want to Continue with Load? YES//<RET>

Will first run the Environment Check Routine, LAMIVTLB

.

Select PackMan function: <RET>

Select MESSAGE Action: IGNORE (in IN basket)// ^<RET>

Select MailMan Menu Option: <RET>

Select Kernel Installation & Distribution System Option: **Installation<RET>**

Select Installation Option: **Install Package(s) <RET>**

Select INSTALL NAME: **LA\*5.2\*12** Loaded from Distribution

4/1/96@07:30: **<RET>**

49 => LA\*5.2\*12 TEST 2

This Distribution was loaded on Apr 01, 1996@07:30:49 with header of  
LA\*5.2\*12

It consisted of the following Install(s): LA\*5.2\*12 **<RET>**

Will first run the Environment Check Routine, LAMIVTLB

Envirnment Check is Ok ---

Install Questions for LA\*5.2\*12

61.38 MICRO INSTRUMENT SET-UP (including data)

Note: You already have the 'MICRO INSTRUMENT SET-UP' File.

Want my data merged with yours? YES// .

61.39 LABORATORY INSTRUMENT CODE (including data)

Note: You already have the 'LABORATORY INSTRUMENT CODE' File.

Want my data merged with yours? YES// .

Want to DISABLE Scheduled Options and Menu Options? YES// NO..

Enter the Device you want to print the Install messages.

You can queue the install by enter a 'Q' at the device prompt.

Enter a '^' to abort the install.

DEVICE: HOME// TELNET

LA\*5.2\*12

-----  
Install Started for LA\*5.2\*12 :

Apr 01, 1996@10:00:05..

Installing Routines:

Apr 01, 1996@10:00:07..

Installing Data Dictionaries: ..

Apr 01, 1996@10:00:09

Installing Data: ..

Apr 01, 1996@10:00:17

Installing PACKAGE COMPONENTS:

## LA\*5.2\*12 Installation Process

Installing PRINT TEMPLATE



Running Post-Install Routine: ^LAMIVTLE

Option [LA VITEK LITERAL] was added to [MICROBIOLOGY] MENU

-----  
-

Option [LA PRINT ORGANISM] was added to [MICROBIOLOGY PRINT] MENU

Option [LA PRINT ANTIBIOTIC] was added to [MICROBIOLOGY PRINT] MENU

Post install completed

Updating Routine file...

Updating KIDS files...

LA\*5.2\*12 Installed.

Install Completed

Post install completed

..

Updating Routine file...

Updating KIDS files...

LA\*5.2\*12 Installed.

Apr 01, 1996@07:31:36..

-----

-

100%		25	50	75	
Complete	+	+			

Install Completed

# Post-Installation Information

## Laboratory Microbiology Staff

### Print DHCP and Vitek Code Entries

1. Print out a list of the DHCP and Vitek antibiotic code entries from the LIC file (#61.39) using the new Antibiotic Code Print [LA PRINT ANTIBIOTIC] option that was added to [MICROBIOLOGY PRINT] MENU by the installation process.

#### **Example:** Antibiotic code entries printout

LABORATORY INSTRUMENT CODE LIST		MAR 25,1996 11:01	PAGE 1
VITEK CODE	VITEK ENTRY	DHCP ENTRY	
-			
amc	Amoxicillin/CA	Amoxicillin	
gm	Gentamicin	GENTMCN	
tax	Cefotaxime	CEFOTAXIME	
mz	Mezlocillin	MEZLOCILLIN	
pip	Piperacillin	PIPERACILLIN	
chl		CHLORAM	
fos	Fosfomycin	Fosfomycin	
FOSFOMYCIN			
dx	Doxycycline	Doxycycline	
DOXYCYCLINE			
e	Erythromycin	ERYTHROMYCIN	
ERYTHROMYCINE			
efx	Enrofloxacin (Vet.)	Enrofloxacin (Vet.)	

2. Print out a list of the DHCP and Vitek organism code entries from the LIC file (#61.39) using the new Organism Code Print [LA PRINT ORGANISM] option that was added to [MICROBIOLOGY PRINT] MENU by the installation process.

#### **Example:** Organism code entries printout

LABORATORY INSTRUMENT CODE LIST		MAR 25,1996 11:10	PAGE 1
VITEK CODE	VITEK ENTRY	DHCP ENTRY	
-			
TODAY	Haemophiluslikebacilli-	HAEMONCHUS, NOS	
gvc	Gramvariablecocci-		
gbv	Gramvariablebacilli-		
gpc	Grampositivecocci-	GRAM POS ROD	
gndip	Gramnegativediplococci-	GRAM NEG ROD	
gnb	Gramnegativebacilli-	GRAM NEG ROD	
gvl		GARBAGE DISPOSAL APP	
ecl	E.colilikebacilli-	ESCHERICHIA COLI	

## Post-Installation Information

bhs	BetahemolyticStrep-	BETA HEMOLYTIC STREP
dl		
camjej	Campylobacterjejuni	CAMPYLOBACTER JEJUNI
dil	diphtheroidlikebacilli	DIPHATHEROIDS

## Editing DHCP Files

The DHCP ENTRY comes from the ETIOLOGY FIELD file (#61.2). The VITEK ENTRY originally comes from the VITEK and is loaded into ETIOLOGY FIELD file (#61.2) by the installation process. If there is a match to the VITEK entry in the ETIOLOGY FIELD file (#61.2) then the LOCAL ORGANISM ENTRY is filled in. A number in the LOCAL ORGANISM ENTRY is not valid; and indicates there was no match or more than one match which could not be resolved. If the organism is one you isolate then decide which entry in the ETIOLOGY FIELD file (#61.2). matches the Vitek entry. If there is no Etiology entry then you must create one and it has to have a SNOMED code. Follow the instructions in the Microbiology Section of the Lab documentation for Etiology File Entries.

**IMPORTANT:** DO NOT change the antibiotic and organism codes on the Vitek or add to the HOST code. This will cause the Literal Interface to become inoperative.

### **LABORATORY INSTRUMENT CODE FILE file (#61.39)**

The LIC file (#61.39) was created to enable editing of the ANTIMICROBIAL SUSCEPTIBILITY file (#62.06) and ETIOLOGY FIELD file (#61.2) for LOCAL ENTRIES.

**Example:** How to edit/add a new antibiotic code using VA FileMan Version 21.0.

Select OPTION:     **ENTER OR EDIT FILE ENTRIES<RET>**

INPUT TO WHAT FILE:     **LABORATORY INSTRUMENT CODE//<RET>**  
 EDIT WHICH FIELD:       ALL//     **ANTIBIOTIC CODE<RET>**             (multiple)<RET>  
       EDIT WHICH ANTIBIOTIC CODE SUB-FIELD:    ALL//<RET>  
 THEN EDIT FIELD: <RET>

Select LABORATORY INSTRUMENT CODE:   **VITEK<RET>**  
 Select ANTIBIOTIC CODE:     tet//<RET>  
       ANTIBIOTIC CODE:     tet//<RET>  
       ANTIBIOTIC INTERPRETATION:   Tetracycline<RET>  
       LOCAL **ANTIBIOTIC** ENTRY:     TETRCLN//

Select LABORATORY INSTRUMENT CODE:   **VITEK<RET>**

```

Select ANTIBIOTIC CODE:   tet//amp <RET>
  Are you adding 'amp' as a new ANTIBIOTIC CODE (the 8th for this LABORATORY
INSTRUMENT CODE)? N<RET>
  (No) ??
Select ANTIBIOTIC CODE: tet// am
  1 am
  2 amc
  3 ams
CHOOSE 1-3: 1<RET>
  ANTIBIOTIC CODE:   am
  ANTIBIOTIC INTERPRETATION:   Ampicillin
  LOCAL ANTIBIOTIC ENTRY:   AMPICLN

```

### Example: How to edit/add a new organism code using VA FileMan Version 21.0.

```

Select OPTION:           ENTER OR EDIT FILE ENTRIES<RET>

INPUT TO WHAT FILE:   LABORATORY INSTRUMENT CODE//<RET>
EDIT WHICH FIELD:     ALL// ORGANISM CODE (multiple)<RET>
  EDIT WHICH ORGANISM CODE SUB-FIELD:     ALL//<RET>
THEN EDIT FIELD:

```

```

Select LABORATORY INSTRUMENT CODE: VITEK<RET>
Select ORGANISM CODE: acibac<RET>
  ORGANISM CODE: acibac//<RET>
  ORGANISM INTERPRETATION:   Acinetobacter baumannii
  Replace
  LOCAL ORGANISM ENTRY:   ACINETOBACTER BAU
  Are you adding 'ACINETOBACTER BAU' as a new ETIOLOGY FIELD (the
4773RD) ???<RET>
Answer with ETIOLOGY FIELD NAME, or SNOMED CODE, or SYNONYM, or
  *BIOCHEMICAL WORKUP
  LOCAL ORGANISM ENTRY:   ACINETOBACTER BAU
  Are you adding 'ACINETOBACTER BAU' as
  a new ETIOLOGY FIELD (the 4773RD)? Y (Yes)
  ETIOLOGY FIELD SNOMED CODE: 7001
  ALREADY AN ENTRY
  ??
  ANSWER MUST BE 2-7 CHARACTERS IN LENGTH
  ETIOLOGY FIELD SNOMED CODE: 7000123

```

```

Select LABORATORY INSTRUMENT CODE:           VITEK<RET>
Select ORGANISM CODE: acibac//e<RET>
  1 e00007
  2 e25922
  3 e35218
  4 e51446
  5 ec
TYPE UP '^' TO STOP, OR
CHOOSE 1-5: 1<RET>
  ORGANISM CODE: e00007//<RET>
  ORGANISM INTERPRETATION: Escherichiacoli//<RET>

```

## Post-Installation Information

LOCAL ORGANISM ENTRY: ESCHERICHIA COLI//<RET>

## LOAD/WORK LIST file (#68.2)

The following example displays how to set up a Load/work List for the Vitek using VA FileMan 21.0 Enter or Edit File Entries option.

### Example

Select OPTION: **ENTER OR EDIT FILE ENTRIES**<RET>

INPUT TO WHAT FILE: **LOAD/WORK LIST**//<RET>

EDIT WHICH FIELD: **ALL**//<RET>

Select LOAD/WORK LIST NAME: `43 VITEK LITERAL<RET>  
 NAME: VITEK LITERAL//<RET>  
 LOAD TRANSFORM: UNIVERSAL//<RET>  
 TYPE: TRAY,CUP//<RET>  
 CUPS PER TRAY: 30//<RET>  
 FULL TRAY'S ONLY: NO//<RET>  
 EXPAND PANELS ON PRINT: <RET>  
 INITIAL SETUP: <RET>  
 VERIFY BY: ACCESSION//<RET>  
 SUPPRESS SEQUENCE #: <RET>  
 INCLUDE UNCOLLECTED ACCESSIONS: NO//<RET>  
 SHORT TEST LIST: <RET>  
 AUTO MICRO EDIT TEMPLATE: ZZTHECULTURESPECIAL//<RET>  
 WKLD METHOD: BACTERIAL ID/SUSC AUTO VITEK//<RET>  
 MAJOR ACCESSION AREA: MICROBIOLOGY//<RET>  
 LAB SUBSECTION: MICROBIOLOGY//<RET>  
 WORK AREA: <RET>  
 DATE OF SETUP: MAR 21,1996//<RET>  
 FIRST TRAY: 1//<RET>  
 STARTING CUP: 1//<RET>  
 LAST TRAY: 1//<RET>  
 LAST CUP: 4//<RET>  
 BUILDING IN PROGRESS: NO//<RET>  
 Select PROFILE: VITEK//<RET>  
 PROFILE: VITEK//<RET>  
 Select TEST: ANAEROBE CULTURE & GRAM STAIN//<RET>  
 TEST: ANAEROBE CULTURE & GRAM STAIN//<RET>  
 SPECIMEN: <RET>  
 BUILD NAME ONLY: NO//<RET>  
 Select TEST: <RET>  
 ACCESSION AREA: MICROBIOLOGY//<RET>  
 Select TRAY #: <RET>  
 Select Specimens to EXCLUDE!: <RET>  
 Select CONTROLS TO BEGIN WORKLIST: <RET>  
 Select CONTROLS TO END WORKLIST: <RET>  
 Select PROFILE: <RET>  
 USER ACCESS AUTHORIZATION: LRVERIFYLABONLY//<RET>  
 Select ADDITIONAL LAB TESTS: <RET>

## AUTO INSTRUMENT file (#62.4)

The AUTO INSTRUMENT file (#62.4) contains all specific information related to each auto instrument in the Laboratory. It is necessary to review and edit this file prior to interfacing the Vitek instrument to run on-line. The following example displays how to add the Vitek Instrument to the AI file (#62.4) using VA FileMan 21.0 Enter or Edit File Entries.

### Example

```
Select OPTION:      ENTER OR EDIT FILE ENTRIES<RET>

INPUT TO WHAT FILE: AUTO INSTRUMENT//<RET>
EDIT WHICH FIELD: ALL//<RET>

Select AUTO INSTRUMENT NAME:      VITEK3<RET>
NAME: VITEK3//<RET>
WKLD METHOD: <RET>
ECHO DEVICE: LSI #1//<RET>
PROGRAM: LAMIVTLP//<RET>
LOAD/WORK LIST: VITEK LITERAL//<RET>
ENTRY for LAGEN ROUTINE: Accession cross-reference           //<RET>
CROSS LINKED BY: +ID//<RET>
*ECHO ALL INPUT: NO//<RET>
METHOD: VITEK//<RET>
DEFAULT ACCESSION AREA: MICROBIOLOGY//<RET>
OVERLAY DATA: YES//<RET>
NEW DATA: D NEW^LASET//<RET>
RESTART: D RESTART^LASET//<RET>
HANDSHAKE RESPONSE: D ^LAMIVTLX//<RET>
ACK TRIGGER VALUE: 5//<RET>
ACK RESPONSE VALUE: 6//<RET>
DIRECT DEVICE: <RET>
Select TEST: ^DEF<RET>
      1  DEFAULT ACCESSION AREA
      2  DEFAULT AUTO MICRO TEST
CHOOSE 1-2: 1<RET>
DEFAULT ACCESSION AREA: MICROBIOLOGY//<RET>
OVERLAY DATA: YES// ^LOAD<RET>
      1  LOAD CHEM TESTS
      2  LOAD/WORK LIST
CHOOSE 1-2: 1 <RET>
LOAD CHEM TESTS: 1//<RET>
Select ALARM TERMINAL: ^<RET>
```



The following example displays how to view the new Vitek entry that was added to AI file (#62.4) using VA FileMan 21.0 Inquire option.

### Example

```

NUMBER: 3                                NAME: VITEK3
ECHO DEVICE: LSI #1                      PROGRAM: LAMIVTLP
LOAD/WORK LIST: VITEK LITERAL            LOAD CHEM TESTS: 1
ENTRY for LAGEN ROUTINE: Accession cross-reference
CROSS LINKED BY: +ID                     *ECHO ALL INPUT: NO
METHOD: VITEK                           DEFAULT ACCESSION AREA: MICROBIOLOGY
OVERLAY DATA: YES                      ACK TRIGGER VALUE: 5
ACK RESPONSE VALUE: 6
DEFAULT AUTO MICRO TEST: CULTURE & SUSCEPTIBILITY
HANDSHAKE RESPONSE: D ^LAMIVTLX         NEW DATA: D NEW^LASET
RESTART: D RESTART^LASET

```

### MICRO INSTRUMENT SET-UP file (#61.38)

The MICRO INSTRUMENT SET-UP file (#61.38) provides another dimension for performing tests on the Vitek. The Vitek only permits one test/accession number. This can be circumvented by entering a number between 0-9 in the prefix field of the LABORATORY TEST file (#60). Employing the logic that the Vitek will only accept up to a 6 digit accession number, and no accession shall exceed 99,999. The prefix number could be placed in front of the 5 digit accession. File (#61.38) also turns on the wild card fields for the Vitek.

**CAUTION:** If you download entries on a patient and that patient already exists in the VITEK database, the "WILD CARD" fields are **OVERWRITTEN** to match the most recent entry. This includes the patient wild card field (pw2) currently being used for physician. If wild cards are used for location, the same thing would happen (overwriting of previous entries for the same patient). There are two concerns: **1)** the epidemiology function in Vitek **CANNOT!** be used because Vitek can enter **ONLY** sort information for **TABLE-DRIVEN** fields and **2)** each time a specimen is downloaded, physician name and patient location for the current specimen overwrites those on previous specimens for the same patient (if they exist). However, the DHCP record would be the official record, but as a back-up, the VITEK database is suspect because the ordering physician (which is required by CAP regulations to be on the report) may not be correct. (Patients could have cultures requested by one physician, get transferred to another location, and have more cultures requested by another physician) For those laboratories wishing to use the epidemiology functions in Vitek, **TABLE-DRIVEN** fields **MUST** be utilized. Therefore, we recommend that the option to use **EITHER** wild card fields **OR** table-driven fields be offered, as well as a mechanism to designate the codes to use for the various downloaded fields during set-up (i.e., FileMan edit by ADPAC).

**NOTE:** Please reference the Bidirectional Computer Interface Operator's Manual, 510710-1, REV 1094 or contact bioMerieux Vitek, 595 Anglum Drive, Hazelwood, MO. 63042-2395, Phone: 314-731-8500, Fax: 314-731-8800 for further instructions.

**NOTE:** If you decide to use the wild card, make certain you change the wild card labels to reflect what you are doing. You will need to change both the data entry (patient data) report to report data.. Be certain to change both the text label and the data label for each wild card field. Delete the labels you will not be using. Follow the instructions given in the bioMerieux or bioLiaison manuals for changing data labels.

The following example displays how to setup Vitek in the MICRO INSTRUMENT SET-UP file (#61.38) using VA FileMan 21.0 Enter or Edit File Entries option.

### Example

Select OPTION: **ENTER OR EDIT FILE ENTRIES<RET>**

INPUT TO WHAT FILE: LABORATORY INSTRUMENT CODE// **MICRO INSTRUMENT SET-UP<RET>**

(1 entry)

EDIT WHICH FIELD: ALL//**<RET>**

Select MICRO INSTRUMENT SET-UP FILE NAME: **VITEK<RET>**

NAME: VITEK//**<RET>**

PREFIX: YES//**<RET>**

Select WILD CARDS ON: YES//**<RET>**

WILD CARDS ON: YES//**<RET>**

w1: LRWRD//?**<RET>**

ENTER ONE

Choose from:

WARD LRWRD

DOC LRDOC

SPEC LRSPEC

TREATING SPECIALTY LRTREAT

BEEPER LRBEEP

w1: LRWRD//**<RET>**

w2: LRDOC//**<RET>**

w3: LRBEEP//**<RET>**

w4: LRSPEC//**<RET>**

w5: LRTREAT//<RET>

## Vitek Literal Verification Process

The following is a representation of the Vitek Literal verification card dependent data as observed with the new DHCP Laboratory Vitek Literal Interface system.

### Example

Select Laboratory DHCP Menu Option: **Microbiology menu<RET>**

```
RB      Results entry (batch)
RE      Results entry
VS      Verification of data by supervisor
VT      Verification of data by tech
WKLD    Review accession workload
        Accessioning, standard (Microbiology)
        Automated Microbiology Menu ...
        Batch accessioning
        Lab statistics menu ...
        Long form accession list for microbiology
        Microbiology print menu ...
        VITEK
        References ...
        Results menu ...
        Short accession list
        Show list of accessions for a patient
        Std/QC/Reps Manual Workload count
        Supervisor menu ...
```

Select Microbiology menu Option: **VITEK**

Vitek Literal verification screen 1

Select auto instrument here: **VITEK JR<RET>**

Select ACCESSION AREA: MICROBIOLOGY//**<RET>**

Accession date: 96//**<RET>** (1996)

Work Load Area: VITEK

Enter the number portion of the Accession: 3161// 3815

There are 7 LAH entries associated with accession 3815

Select one of the following:

```
1      OK
2      NO
```

Enter response: OK//**<RET>**

ACC # (3815) FEB 15,1996@10:57:48

DEMO,DENNIS S SSN: 000-00-0114 LOC: LAB  
Specimen: FECES Sample: STOOL

GRAM STAIN 4+ nr  
3 + nr  
2+ gram negative rod

Is this the correct patient/specimen? ? Yes//<RET> (Yes)

DEMO,DENNIS S 000-00-0114  
MI 96 3815

FECES STOOL

- 1 GRAM STAIN
- 2 CULTURE & SUSCEPTIBILITY
- 3 CAMPYLOBACTER CULTURE

Please enter the test number(s) or : ALL//<RET>

ENTER QUANTITY FOR ( KLEBSIELLA PNEUMONIAE ) : 4+ // (4+)

4+ ? Yes// (Yes)

COMMENT: //

ENTER QUANTITY FOR ( ESCHERICHIA COLI ) : 4+ // (4+)

4+ ? Yes// (Yes)

COMMENT: //

Isolate (1 )

KLEBSIELLA PNEUMONIAE

CARD gns-f3

AMIKACN	<=2	S	A
CEFAZOLIN	<=8	S	A
CEFOPERAZONE	<=8	S	R
CEFOTAXIME	<=4	S	A
CEFOXITIN	8	S	A
CEFTAZIDIME	<=8	S	A
CEFUROXIME	<=4	S	R
CEPHALOTHIN	4	S	A
NITROFURANTOIN	<=32	S	A
TOBRMCN	<=0.5	S	A

Enter RETURN to continue or '^' to exit:

Isolate (1 )

KLEBSIELLA PNEUMONIAE

CARD gns-f6

AMPICILLIN	>=32	R	A
AMPICILLIN/SULBACTAM	16	I	A

## Post-Installation Information

AZTREONAM	<=8	S	R
CIPROFLOXACIN	<=0.5	S	A
IMIPENUM	<=4	S	R
MEZLOCILLIN	<=16	S	R
OFLOXACIN	<=1	S	A
PIPERACILLIN	<=8	S	A
TICARCILLIN	>=256	R	R
TICARCILLIN/CA	<=16	S	R
TRMSULF	<=10	S	A

Enter RETURN to continue or '^' to exit:

Isolate (2 )

ESCHERICHIA COLI

CARD gni

Enter RETURN to continue or '^' to exit:

Isolate (2 )

ESCHERICHIA COLI

CARD gns-f3

AMIKACN	<=2	S	A
CEFAZOLIN	<=8	S	A
CEFOPERAZONE	<=8	S	R
CEFOTAXIME	<=4	S	A
CEFOXITIN	<=2	S	A
CEFTAZIDIME	<=8	S	A
CEFUROXIME	<=4	S	R
CEPHALOTHIN	8	S	A
NITROFURANTOIN	<=32	S	A
TOBRMCN	<=0.5	S	A

Enter RETURN to continue or '^' to exit:

Isolate (2 )

ESCHERICHIA COLI

CARD gns-f6

AMPICILLIN	>=32	R	A
AMPICILLIN/SULBACTAM	16	I	A
AZTREONAM	<=8	S	R
CIPROFLOXACIN	<=0.5	S	A
IMIPENUM	<=4	S	R
MEZLOCILLIN	>=256	R	R
OFLOXACIN	<=1	S	A
PIPERACILLIN	>=256	R	A
TICARCILLIN	>=256	R	R
TICARCILLIN/CA	<=16	S	R
TRMSULF	<=10	S	A

Enter RETURN to continue or '^' to exit:

BACT RPT STATUS: ??

Choose from:

F FINAL REPORT

P PRELIMINARY REPORT

BACT RPT STATUS: P PRELIMINARY REPORT

Select GRAM STAIN: 2+ gram negative rod //  
 Select BACT RPT REMARK:

DEMO,DENNIS S SSN: 000-00-0114  
 MI 96 3815 STOOL FECES

1. KLEBSIELLA PNEUMONIAE (4+ )		2. ESCHERICHIA COLI (4+ )		
:	:	:	:	
SUSC	INTP	SUSC	INTP	
AMIKACN	<=2 S	<=2 S		\$\$\$\$
AMPICILLIN/SUL	16 I	16 I		\$
AMPICILLIN	>=32 R	>=32 R		\$
AZTREONAM	<=8* S*	<=8* S*		\$\$\$\$
CEFAZOLIN	<=8 S	<=8 S		\$
CEFOPERAZONE	<=8* S*	<=8* S*		\$
CEFOTAXIME	<=4 S	<=4 S		\$
CEFOXITIN	8 S	<=2 S		\$
CEFTAZIDIME	<=8 S	<=8 S		\$
CEFUROXIME	<=4* S*	<=4* S*		\$
CEPHALOTHIN	4 S	8 S		\$
IMIPENUM	<=4* S*	<=4* S*		\$
MEZLOCILLIN	<=16* S*	>=256*R*		\$
NITROFURANTOIN	<=32 S	<=32 S		\$
OFLOXACIN	<=1 S	<=1 S		\$

PRESS RETURN FOR MORE

DEMO,DENNIS S SSN: 000-00-0114  
 MI 96 3815 STOOL FECES

1. KLEBSIELLA PNEUMONIAE (4+ )		2. ESCHERICHIA COLI (4+ )		
:	:	:	:	
SUSC	INTP	SUSC	INTP	
PIPERACILLIN	<=8 S	>=256 R		\$\$\$\$
TICARCILLIN	>=256*R*	>=256*R*		\$\$\$\$
TICARCILLIN/CA	<=16* S*	<=16* S*		\$\$\$\$
TRMSULF	<=10 S	<=10 S		\$

('E'dit data, 'C'omments, 'O'rganism 'W'orklist) // ^

REMOVING ^LR DATA







# Troubleshooting

1. **Problem:** No Data in ^LA

*Possible Solutions:* Check Vitek protocol setting  
Check wiring connections

2. **Problem:** No Data in ^LAH

*Solutions:* Loadlist not defined  
Loadlist not defined in LIC file (61.39)  
Organism defined in LIC file (61.39)  
Antibiotic defined in LIC file (61.39)

3. **Problem:** Susceptibility not showing up on reports

*Solution:* Check the ANTIMICROBIAL SUSCEPTIBILITY file (#62.06) to see if the MIC from the Vitek matches that in the Vitek file (i.e. Vitek=0.5, file=.5, file must be changed to 0.5).